

# CU-120

100% CONTENT UNIFORMITY


PHARMA **TECHNOLOGY** 



API's



WEIGHT

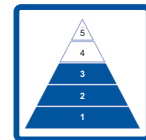

≤ OEB3

New and innovative in-line solution designed to ensure the quality of each individual tablet or capsule, by combining two main control methods: API's fraction and weight.

The **CU-120** analyses individual dosage from each single tablet or capsule up to a speed of 120.000 tablets per hour and rejects all non-compliant products.

The ideal tool for Content Uniformity Inspection and RTD (Real Time Decision).

---

**ADVANCE BY TECHNOLOGY**


# 3D MICROWAVE RESONATOR FOR WEIGHT MEASUREMENT

Each product is inspected by the multipoint **NIR-SRS** “near-Infrared spatially resolved technology” **probe to predict APIs concentration**, and a **3D microwave resonator** allowing to **measure individual weight**. Various extensive repeatability **tests campaigns and industrial case studies have shown that the predictions are within the predicted error boundaries**.

The unit include an **encapsulated weighcell**, or a **highly accurate 4-P tester** (measuring weight, thickness, hardness, and diameter), for automatic calibration of the measurement instrument. A linear regression is performed to warranty that the accuracy remains inside the **machine accuracy of +/-1mg (3 Sigma)**.

Data are processed and analysed by hi-speed processors and stored in a **user-friendly “step by step” CFR21 acquisition software**. Data can easily be **exported to an USB memory device**, and **process data can easily be exported to an up-stream machine or software through OPC UA** (e.g. connection to a SIPAT infrastructure or a third-party process equipment).

NIR-SRS probe and 3D microwave resonator

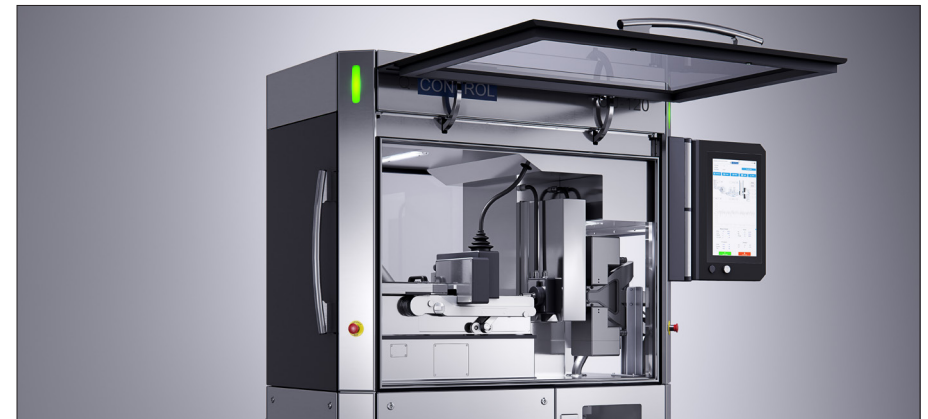

Embedded 4-P tester



**WEIGHT**




**THICKNESS**



**HARDNESS**



**DIAMETER**

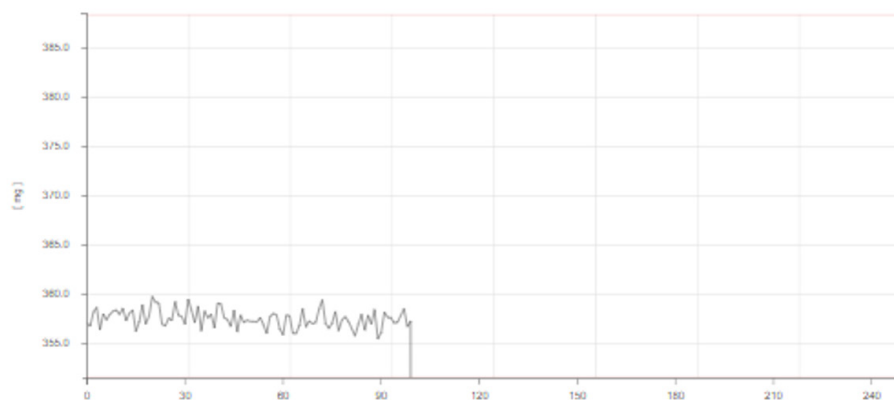

## Measured weight [mg]

Actual	mg	357.27
Count	#	100
Sampled	#	0
Out of spec.	#	0

## Range

Min	mg	355.49
Max	mg	359.78
Average	mg	357.53

## Throughput

Actual	#/sec	0
Average	#/sec	0
Total	#/sec	0

## Deviation

Sabs	mg	0.89
Srel	%	0.25

# NIR SRS TECHNOLOGY SPATIALLY RESOLVED SPECTROSCOPY FOR API'S FRACTION PREDICTION

Equipped with the fasted NIR-SRS technology and an highly precise 3-D microwave weight measurement station this equipment is geared to provide you a **very precise Content Uniformity prediction**, and to remove 100% of non-compliant products from good products, **allowing RTD (Real Time Decision)**

The machine can also be used for **R&D developments, batch recovery/saving, mix-up and counterfeit products.**



The unit has been developed with an utmost care to the mechanical design allowing **fast changeover and affordable format changes.**

An utmost care has been taken to **simplify the equipment to allow use in a production environment.**

Designed to comply with the European Union's CE Guidelines, they are supplied with a manual, safety instructions and wiring diagram.

**Our company can assist you for installation on site and help you to develop chemometric models.**

- 100% Content Uniformity test of your products
- Permit Real Time Decision
- Ideal for proving final quality on batch or CM
- Shorten your reject strategy process
- Test physical attributes of your tablets based on conventional sampling methods (weight, thickness, hardness, and length/diameter)
- Inspection machine for batch recovery, prove global batch quality, detect counterfeit product

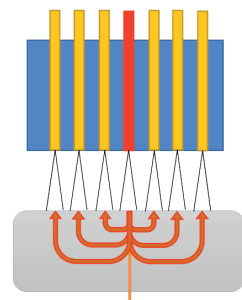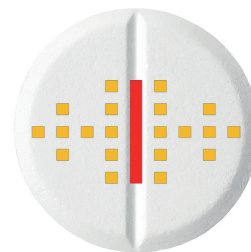


- Light Source
- Collection channels

NIR-SRS multipoint probe



# ADVANCE BY TECHNOLOGY

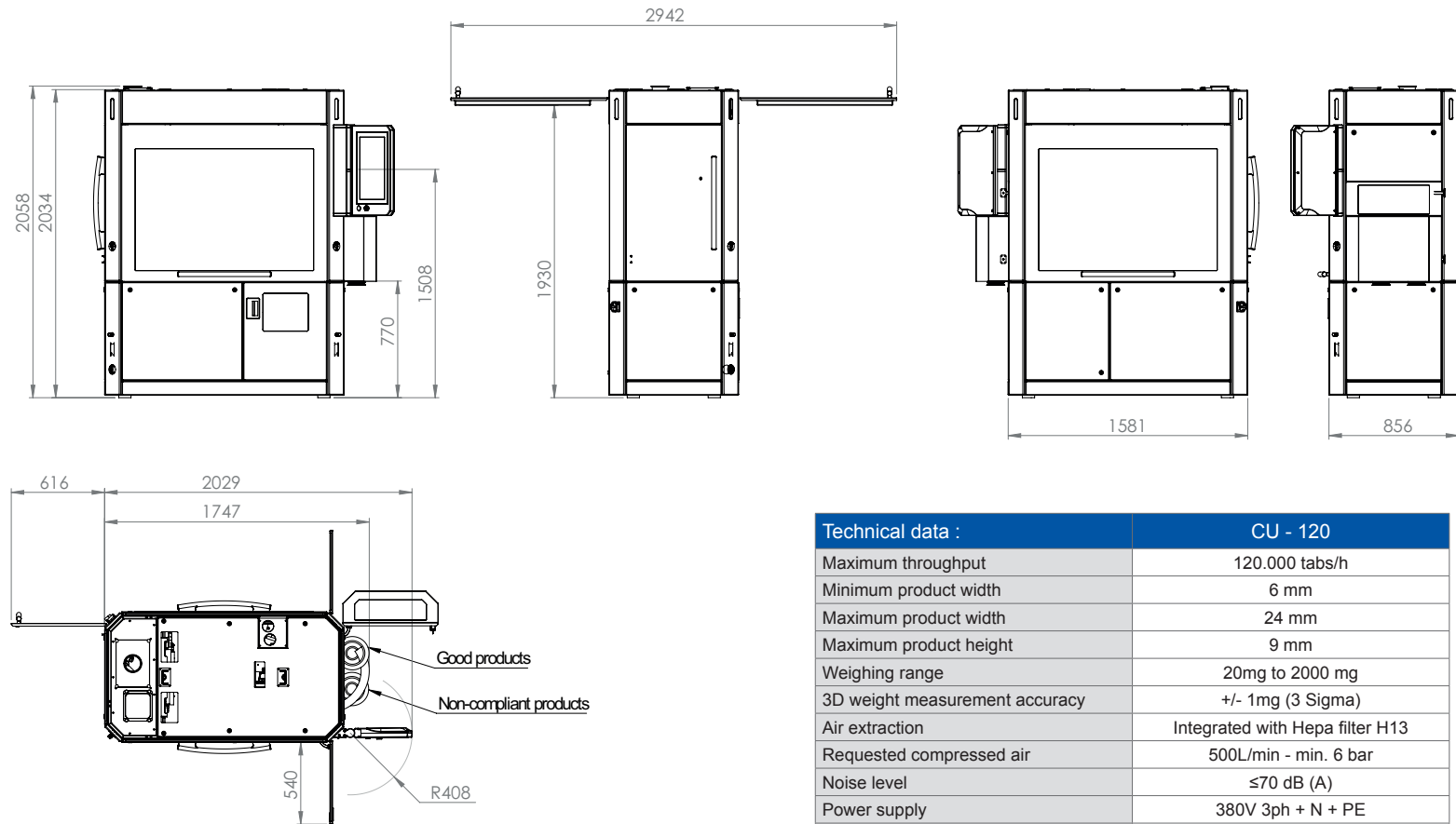

Technical data :	CU - 120
Maximum throughput	120.000 tabs/h
Minimum product width	6 mm
Maximum product width	24 mm
Maximum product height	9 mm
Weighing range	20mg to 2000 mg
3D weight measurement accuracy	+/- 1mg (3 Sigma)
Air extraction	Integrated with Hepa filter H13
Requested compressed air	500L/min - min. 6 bar
Noise level	≤70 dB (A)
Power supply	380V 3ph + N + PE
Frequency	50 Hz / 60 Hz
Polymer contact parts	FDA approved

PHARMA **TECHNOLOGY** 

Rue Graham Bell 8  
 B-1402 Thines (Nivelles) - Belgium  
 Phone : +32 (0)67 70 13 00  
 E-Mail : info@pharmatec.be